



## Microtest – Main window minimizes at startup

This problem can be caused because one of the files that stores various parameters about the system has become corrupted. Fortunately there is a simple way to resolve this which is described below:

1. Open “My Computer”
2. Navigate to “C:\ProgramData\deben”
  - Note: the ProgramData folder can sometimes be hidden. To show this folder:
    - Press “Alt-T” within Windows explorer to open the Tools menu
    - Click “Folder options”
    - Click on the “View tab”
    - Tick the option that says “Show hidden files, folders and drives”

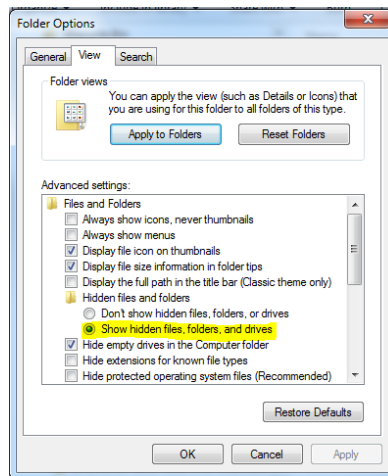

3. In the “Deben” folder you should see two files: mt\_tt.par and winpos.deb. Delete the winpos.deb file.

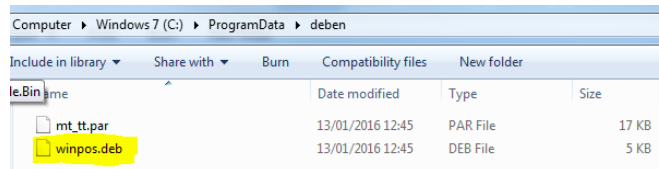

4. Restart the Microtest software. The main window should now appear normally.

If this problem happens regularly, you can permanently disable this function by going into the Module Setup window, clicking on the Global tab and unticking the “Remember Window Positions” checkbox. This should mean that the “winpos.deb” file is no longer created or referenced.

